

Human Anti-PEG IgG ELISA kit (MAK-0019)

This product is for research use only and is not intended for diagnostic use.

PRODUCT INFORMATION

Cat No. MAK-0019

Storage Indirect ELISA

Principle This kit is for research use only. Under no circumstances may it be used for diagnostic purposes.

Intended Use Attachment of polyethylene glycol (PEG) chains to therapeutic biologic agents, a process referred to as PEGylation, prolongs the circulating half-life of the modified protein by slowing proteolytic degradation and by masking it from the immune system. However, it has been reported that repeated injections of PEGylated proteins can induce anti-PEG antibodies that increase the rate of clearance and decrease drug efficacy (accelerated blood clearance, or ABC phenomenon). To aid research in this key area, we have developed a human anti-PEG IgG ELISA kit.

Introduction The assay uses immobilized mono-mPEGylated BSA (20 kDa PEG chain) as the capture antigen (coated on microtiter wells) and horseradish peroxidase (HRP) conjugated anti-human/monkey IgG monoclonal antibody (LDI clone 6D11-21) for detection. Serum and plasma samples are diluted and incubated alongside standards in the microtiter wells for 1 hour. The wells are subsequently washed. HRP conjugate is added and incubated for 45 minutes. Anti-PEG IgG molecules are thus sandwiched between immobilized PEG and the detection antibody conjugate. The wells are then washed to remove unbound HRP-labeled antibodies. TMB reagent is added and incubated for 20 minutes at room temperature. This results in the development of a blue color. Color development is stopped by the addition of Stop Solution, changing the color to yellow. Optical density is measured spectrophotometrically at 450 nm. The concentration of anti-PEG IgG is

proportional to the absorbance at 450 nm and is derived from a standard curve.

Principle of the assay It has been reported that repeated injections of PEGylated proteins can induce anti-PEG antibodies that increase the rate of clearance and decrease drug efficacy (accelerated blood clearance, or ABC phenomenon). To aid research in this important area, we have developed a anti-PEG ELISA kit. Used for new drug research and development

Storage Note The reference stock, HRP conjugate and the PEG-BSA coated plate should be stored at -20°C. All remaining kit components should be stored at 4°C. The microtiter plate should be kept in a sealed bag with desiccant. Kits will remain stable for six months from the date of purchase provided that the components are stored as described.

Materials MATERIALS
Materials provided with the kit:

- PEG-BSA coated plate (12 x 8-wells)
- Anti-IgG HRP Stock (lyophilized)
- Reference Stock (lyophilized)
- 20x HRP PEG Wash: PEGW50-20, 50 ml
- HRP PEG Diluent: PEGD50-1, 50 ml
- TMB: TMB11-1, 11 ml
- Stop Solution: SS11-1, 11 ml

General instructions Please read and understand the instructions thoroughly before using the kit. All reagents should be allowed to reach room temperature (25°C) before use. The wash procedure is critical. Insufficient washing will result in poor precision and falsely elevated absorbance readings. Use only the wash solution and dilution buffer provided with the kit. PEG and PEGylated compounds are found in buffers conventionally used in ELISA's and cannot be used with this kit. Kits are validated using plate shakers set at 150 rpm and 25°C. Performance of the assay at lower temperatures and/or mixing speeds will result in lower absorbance values. Optimal results are achieved if at each step, reagents are pipetted into the wells of the microtiter plate within 5 minutes.

Diluent The diluent (PEGD50-1) is formulated for measurement PEG antibodies. It is supplied ready to use. DO NOT substitute other buffers.

Sample Preparation	In studies at Mabioway Co., Ltd., we found that anti-PEG IgG levels in human samples ranged from undetectable to 3750 u/ml. Optimal dilutions must be determined empirically. However, we suggest that samples initially be diluted 20-fold. To avoid matrix effects, do not test dilutions less than 20-fold (i.e., 10-fold).
HRP Anti-PEG Conjugate	Approximately 15 minutes before needed, reconstitute the lyophilized HRP conjugate as directed on the vial label and mix gently. Dilute as described on the vial label to give the working conjugate solution. The reconstituted conjugate stock must be stored at -20°C in a sealed vial if future use is intended.
Wash Solution	The wash solution is provided as a 20x stock. Prior to use, dilute the contents of the bottle (50 ml) with 950 ml of distilled or deionized water.
Standards	<ol style="list-style-type: none"> 1.The anti-PEG IgG standard is provided as a lyophilized stock. Reconstitute as described on the vial label. 2.Label 8 polypropylene or glass tubes as 100, 50, 25, 12.5, 6.25, 3.13, 1.56 and 0.78 u/ml². 3.In the tube labeled 100 u/ml prepare the 100 u/ml standard as detailed on the stock vial label. 4.Dispense 250 ul of diluent into the remaining tubes. 5.Prepare a 50 u/ml standard by diluting and mixing 250 ul of the 100 u/ml standard with 250 ul of diluent in the tube labeled 50 u/ml. 6.Similarly prepare the remaining standards by serial dilution.
Procedure	<ol style="list-style-type: none"> 1.Secure the desired number of coated wells in the holder. 2.Dispense 100 ul of standards and diluted samples into the wells (we recommend testing in duplicate). 3.Incubate on a plate shaker at 150 rpm/25°C for 1 hour. 4.Aspirate the contents of the microtiter wells and wash the wells five times with 1x wash solution using a plate washer (400 ul/well). 5.Strike the wells sharply onto absorbent paper to remove all residual wash solution. 6.Add 100 ul of diluted HRP conjugate into each well. 7.Incubate on a plate shaker at 150 rpm/25°C for 45-minutes. 8.Wash as detailed above. 9.Dispense 100 ul of TMB into each well. 10. Incubate on a plate shaker at 150 rpm/25°C for 20 minutes. 11. Stop the reaction by adding 100 ul of stop solution to each well.

12. Gently mix. It is important to make sure that all the blue color changes to yellow.
 13. Read the optical density at 450 nm with a microtiter plate reader within five minutes³.
-

Important Tips

/

Results

1. Using curve fitting software, construct a standard curve by plotting absorbance values of the standards versus \log_{10} of the concentration.
 2. Fit the standard curve to a four-parameter logistic regression (4PL) equation (x axis = \log_{10} concentration) and determine the concentration of the samples from the standard curve (remember to derive the antilog).
 3. Multiply the derived concentration by the dilution factor to determine the actual concentration in the samples.
 4. If the A_{450} values of samples fall outside the standard curve, samples should be diluted appropriately and re-tested.
-